

IMPORTANT INFORMATION TO READ and RETURN

Installation Requirements for a Masterclave 09 Media Preparator

Thank you for choosing one of our products for your laboratory. To enable our engineers to perform an efficient, trouble-free installation please study, complete and fax this form to us on **01274 531197**. Should you have any questions, please do not hesitate to contact us, as we are here to help. When we have received the completed form, our Service Department will contact you to arrange a mutually convenient installation date.

The following information represents the ideal requirement.
Please contact us IMMEDIATELY if your intended location does not match this specification.



Access Requirements

The unit is designed to fit through a door opening of 700mm (27.5") but check size of lifts, stairs etc.

☐

Space Requirements

The weight of the equipment is 60kg. The bench allocated must be flat, level and of sufficient size to support the base fully.

The unit should be sited as close as possible (preferably to the right) of the pourer stacker

External Dimensions

Depth	Width	Height
mm	mm	mm
435	500	590

☐

Mains Requirements

Electricity Supply

230V +/- 10% AC, single phase

Socket

1 x Three pin, 13 amp

☐

Water

Maximum pressure – 2 bar (25psi) Flow rate – 4 to 6 litres per minute.

Maximum temperature - 20°C

☐

Connection

3/4 " bsp male (washer connection) fitting or a 1/2 " (15mm) barb. In both cases, an isolation tap is recommended.

☐

Drain

1 3/4 " (30mm) or 1 1/2 " (38mm) waste trap or a sufficiently large bore waste pipe to accommodate two 1" (25mm) hoses.

☐

Validation is highly recommended for each load to be utilised in order to:

- ensure that the required temperatures and times throughout the load are achieved
and
- determine cycle settings.

As Don Whitley Scientific is **UKAS accredited**, we can offer you validation and calibration of your heat sterilisation equipment.

If you would like to discuss this, please call us on 01274 595728.

Delivery and installation are free of charge (unless otherwise agreed). If our engineers are unable to install the unit and a return journey is necessary, **a charge may be made**.

It is essential that this form is completed and returned to avoid delay to your installation.

THANK YOU FOR THINKING WHITLEY

Signature

Title

Print Name

Establishment